



Shipping Refrigerated Embryos



- Completely fill 0.5ml microfuge tube with 0.6ml M2 medium at room temperature.
- Load 30-40 embryos into each microfuge tube, dispensing the embryos towards the bottom of the tube.

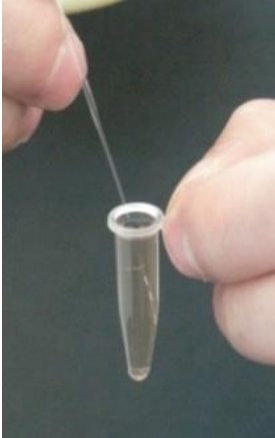


- Seal tube with parafilm.
- Place the tube containing the embryos into biotube which is supplied within the cold transportation kit.

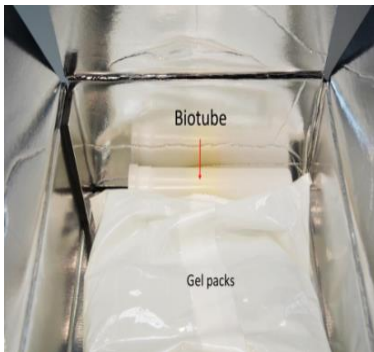

- Place biotube into aluminium lined box (room temperature).
- Place two gel cool packs (room temperature) into aluminium box, surrounding biotube.



- Seal aluminium box with sellotape.
- Place aluminium box into polystyrene container following assembly instructions.

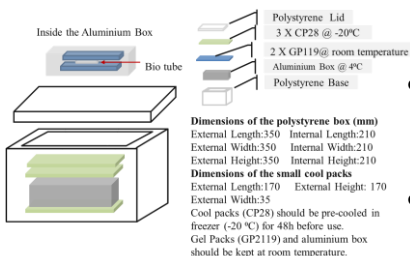

- Seal polystyrene container with packing tape.
- Place polystyrene container into cardboard box.
- Send package via regular delivery service.

